



Department of Chemistry
NMR Facilities
Director: Dr. Carlos A. Steren

The DEPTQ EXPERIMENT

DEPTQ is similar to the DEPT experiment but it also allow us to detect non protonated carbons. Below, the different carbon signals observed on a DEPTQ and on a DEPT are detailed. Blue and red indicate signal phase up or down respectively.

DEPTQ 45°
C, CH, CH₂, CH₃

DEPT 45°
CH, CH₂, CH₃

DEPTQ 90°
C, CH

DEPT 90°
CH

DEPTQ 135°
C, CH, CH₂, CH₃

DEPT 135°
CH, CH₂, CH₃


Setting up the DEPTQ experiment.

DEPTQ is available on the Varian 500 and Varian 600 spectrometers.

1.- Call the “HYDROGEN” experiment and then call the “CARBON” experiment. Both experiments in the same experiment page. (This is necessary for both DEPT and APT due to a bug in VNMRJ.)


2.- Call the DEPT experiment.

If DEPT or APT are called directly (skipping step 1), an alert sign appears at the bottom of the VNMRJ window. This indicates that the experiment was not set up properly (see the figure below).

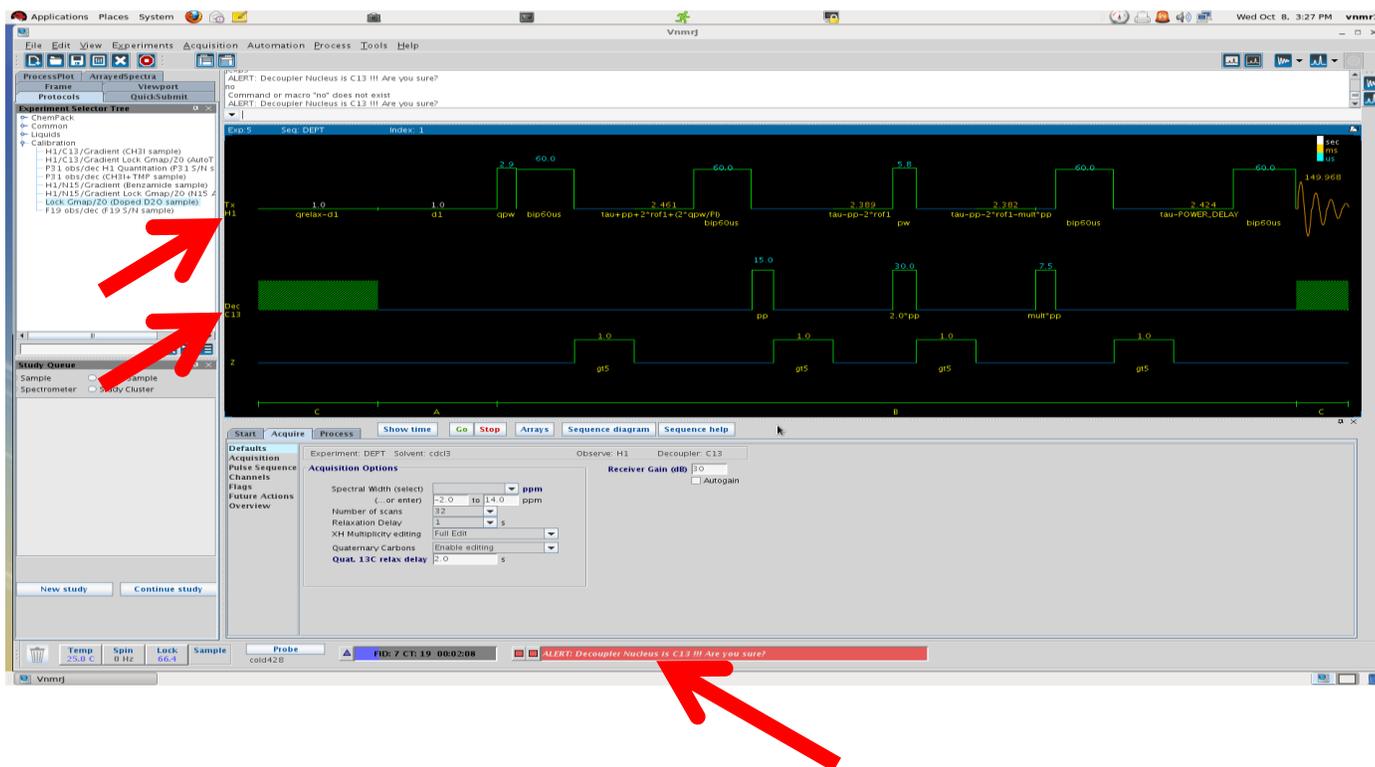


Also, the pulse sequence shows that the experiment will detect H1 and that the decoupler channel is set to ^{13}C . This is wrong!

When the “CARBON” sequence is loaded before the DEPT, there is no alert sign and, the DEPT pulse sequence detects ^{13}C and it decouples ^1H , as shown in the figure below.

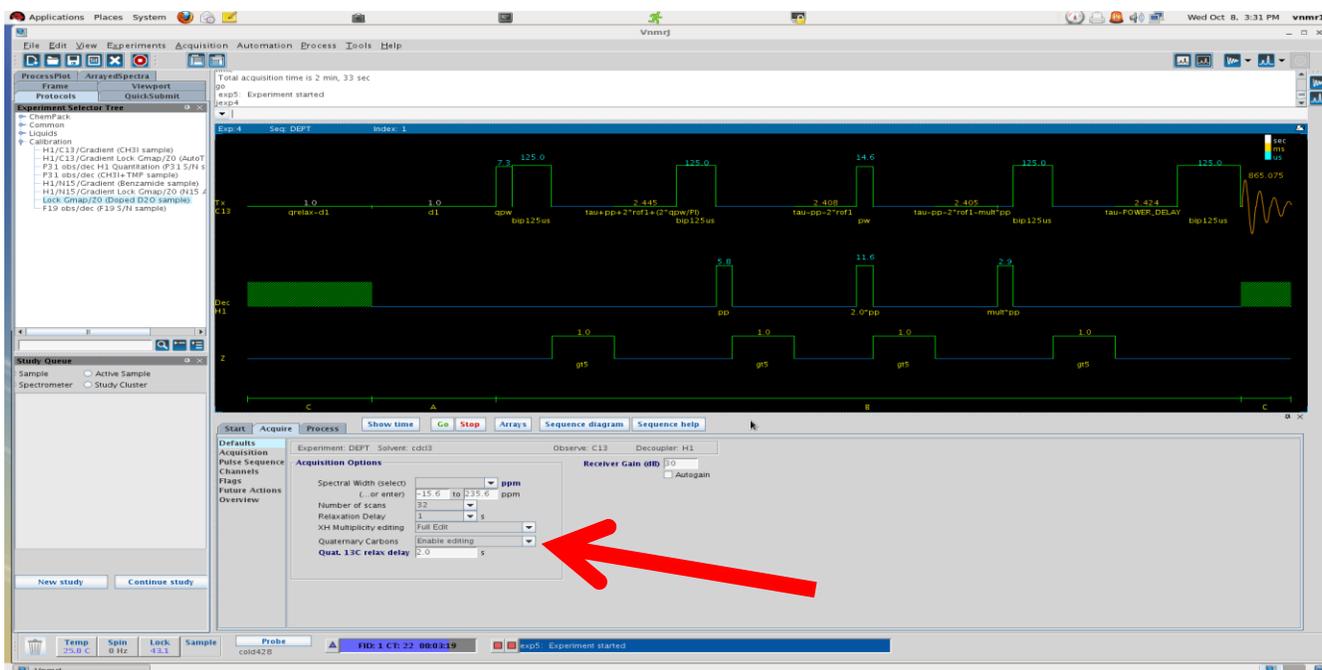

DEPTQ and DEPT options

The default DEPT is now the DEPTQ. To set up the old DEPT, on the “Quaternary Carbons” line, select the option “Suppress”, as shown in the next figure (red arrow).

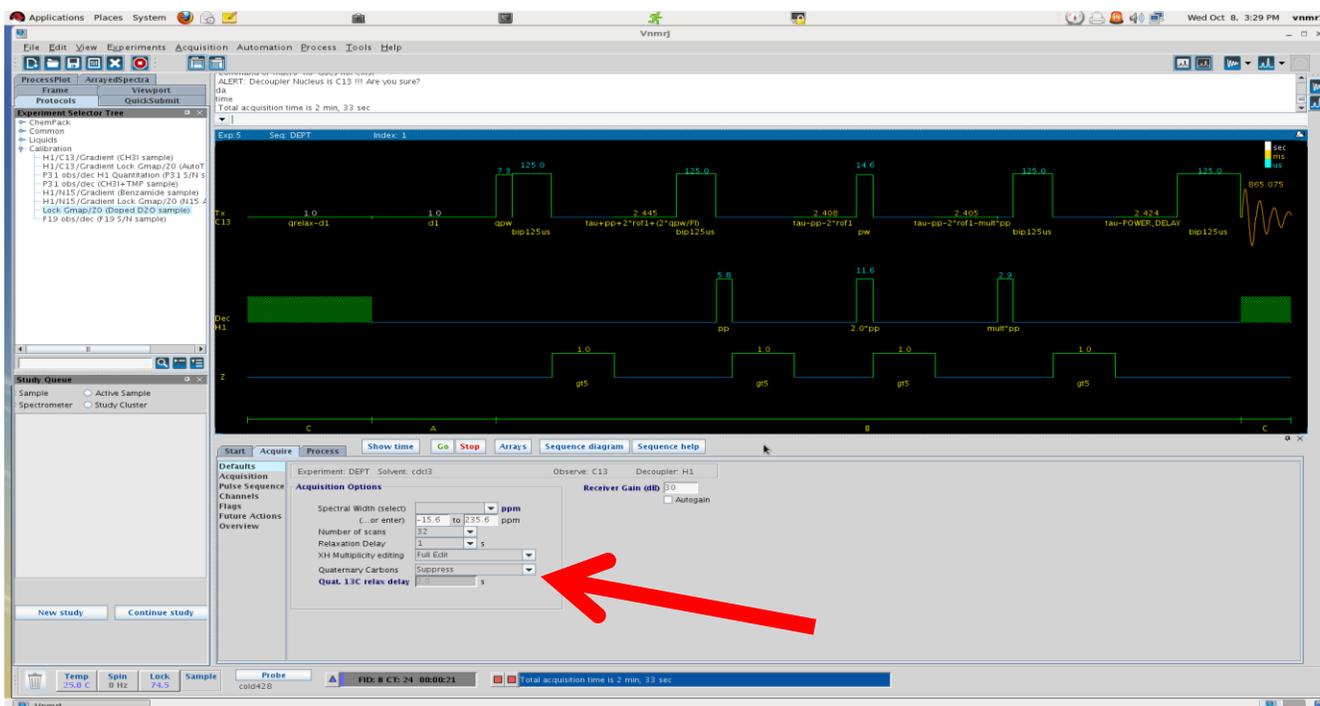

To run only a 90° spectrum, on the “XH multiplicity” line select the “XH only” option. For a 135° spectrum select the “XH/XH3 up & XH2 down” option.

Full editing of carbon spectra:

The DEPTQ generates 8 spectra while the DEPT generates 4 spectra. These spectra can be processed with MestreNova to yield “full carbon edited spectra”.

1.- Open the DEPTQ or DEPT spectra on MestreNova. The software processes the DEPTQ automatically and displays the “full carbon edited spectra” shown below.

